



## **FE-340-H1-PAR-45 - Operating manual**

### **1. Turning the unit ON**

- Press the **ON** button to turn the unit ON – “ON” will appear on display.

### **2. Turning the unit OFF**

- Press the **OFF** button to turn the unit OFF – “OFF” will appear on display.

### **3. Setting temperature**

- Press the **(+)** or **(-)** buttons – the set-point temperature will appear on display.
- Adjust the set-point temperature using the **(+)** or **(-)** buttons.

---

### **4. Technician settings – offset for temperature readings calibration and protective high limit for heating.**

- Press and hold the empty button (without text) for 20 seconds until the offset - a number between -6...+6 appears on display.
- Adjust the offset temperature using the **(+)** or **(-)** buttons. The offset is used for calibration of the measured temperature when needed.
- Press the empty button again – 44, 45 or 46 will appear on display.
- Using the **(+)** and **(-)** buttons adjust the protective high limit for heating. When the temperature reaches the limit, the heat element will stop.
- Press the empty button again to return to normal display.